



3500 Winton Place
Rochester, NY 14623
585-442-0930 (phone)
585-442-0934 (fax)
www.litronlabs.com

Mouse MicroFlow[®] Micronucleus Analysis Kit Biological Standard Set-up Data

The mouse biological standards (Malaria Biostandard, Negative Control, Positive Control, and CD71-Negative) supplied in this MicroFlow[®] Kit, have been carefully prepared to assist you in optimally configuring your flow cytometer for micronucleus scoring.

Upon receipt of the biological standards, immediately place them upright in a -75 °C to -85 °C freezer. Each of the enclosed vials contains approximately 1.0 ml.

The Negative Control and Positive Control biological standards have been analyzed at Litron Laboratories. Your measurements should approximate the values below:

<u>Biological Standard</u>	<u>% RETs</u>	<u>% MN-RETs</u>
Negative Control Lot # MN1012	1.41 to 2.20	0.12 to 0.27
Positive Control Lot # MP1012	0.77 to 1.20	1.42 to 3.19

If you have any questions regarding these procedures or those described in the manual, please contact Litron Laboratories at 585-442-0930 or by email to info@litronlabs.com.